

OPERATING MODE RANGES

- d:** multidispensing, TipGuide makes the best tip choice on behalf of the user
- dST:** multidispensing, user selects the tip size (see table below!)

KEYBOARD

Select



- Press to go to the Main Modes
- Press to go to the previous step
- Activates the speed function

Arrow up/down



- Press to select the mode or make setting changes (volume, aliquots, speed, etc.)

Enter



- Press to confirm the settings

OPERATION BUTTONS

Start button



- Triggers the aspiration and dispensing operations
- Confirms the selected tip

Tip ejection buttons



- Press to remove tip

Reset button



- Press to reset the pipettor

TIP SELECTION WITH INCREMENTS (for dST mode)

Tip Cat. No.	Tip volume (ml)	Volume min/max (µl)	Increment (µl)	Number of dispensing Min/max
40007-0000	0,1	d: 1/100	d: 0,2	1/100
40007-0001	0,2	d: 2/200	d: 0,4	1/100
40007-0003	0,5	d: 5/500	d: 1	1/100
40007-0004	1,0	d: 10/1000	d: 2	1/100
40007-0005	2,5	d: 25/2500	d: 5	1/100
40007-0006	5,0	d: 50/5000	d: 10	1/100
40007-0007	10,0	d: 100/10 000	d: 20	1/100
40007-0009	25,0	d: 500/25 000	d: 50	1/50
40007-0010	50,0	d: 1000/50 000	d: 100	1/50

PROGRAMMING MAP

